

# **INDUSTRIAL HYGIENE REPORT**

## **Control of Formaldehyde**

### **In the Workplace**

**May 2011**

**For:** Sentry Air Systems, Inc.  
Attn.: Omar Ilsever, Sales Manager  
6999 W. Little York, Ste. P1  
Houston, TX 77040

**Prepared by:** Robert F. Adams  
RF ADAMS & ASSOCIATES, INC.  
12111 Millstream Way  
Houston, TX 77041

**Date of tests:** April 08 and April 27, 2011

## Control of Formaldehyde in the Workplace

**Introduction:** Formaldehyde is a gas that is sometimes found in emissions from particle board or from other products made from formaldehyde such as urea-formaldehyde or phenol-formaldehyde polymers. It has been reported<sup>1</sup> that some consumer products used to straighten or treat human hair may contain or release formaldehyde during use. Formaldehyde has a penetrating, irritating odor. It is highly reactive and readily polymerizes to form a trimer (trioxymethylene) or the solid, paraformaldehyde. Formaldehyde is often marketed as a 37% (by weight) water solution called formalin. It is usually stabilized by the addition of 8-10% methanol. Formaldehyde is a confirmed carcinogen and workers' exposure is regulated by OSHA. The eight-hour TWA exposure should not exceed 0.75ppm and the 15-minute short-term exposure limit (STEL) has been set at 2ppm.

**Purpose:** The purpose of this experiment was to demonstrate the effectiveness of the Sentry Air Systems Model 300 *Winged Sentry™* with Lid (WSL) in collecting and removing formaldehyde gas produced by the evaporation of a measured amount of liquid formalin. The Model 300 WSL was equipped with an activated charcoal filter specially treated for the removal of aldehydes for one test and a standard activated charcoal filter for another test. There are two WSL models available (Model SS-200 WSL and Model SS-300 WSL); the latter was used in this project.

**Discussion:** Formaldehyde is best captured near its source and the Model 300 WSL is very efficient in capturing and removing formaldehyde before it mixes with the surrounding workplace air. Formaldehyde that escapes into the room air can be a health concern for personnel since it has poor warning properties and air concentrations could exceed the recommended limit before its odor is detected. For purposes of this experiment a measured amount of liquid formalin was distributed drop-wise from a burette onto several cotton balls placed in a shallow pan. The cotton provided a large surface area to facilitate the evaporation of the formaldehyde solution. When the Model 300 WSL was turned on, a stream of air was drawn across the cotton balls and into the Sentry fume extractor.

<sup>1</sup>[http://articles.chicagotribune.com/2011-04-12/kkmk/chi-hair-straightening-products-could-release-formaldehyde-20110412\\_1\\_brazilian-blowout-oregon-s-osh-formaldehyde](http://articles.chicagotribune.com/2011-04-12/kkmk/chi-hair-straightening-products-could-release-formaldehyde-20110412_1_brazilian-blowout-oregon-s-osh-formaldehyde)

**Experimental:** On April 8, 2011, an activated carbon filter treated for the removal of aldehydes was prepared and installed in the Model 300 WSL. The air flow rate with this specific filter was measured to be approximately 250 cfm (cubic feet per minute). 4.0 ml of formalin (containing 1.48 grams of formaldehyde) was dispensed from a 50-ml burette and evaporated into the air stream over a 30-minute period. Air samples were collected during the experiment using SKC 226-119 sorbent tubes and battery-powered sampling pumps. The pumps were set to pull about 100 cc/min. and were pre- and post-calibrated. All personnel involved in the sampling project wore half-face, air-purifying respirators equipped with an organic vapor cartridge while formalin was being handled. The operator also wore eye protection. Samples were taken at the inlet and outlet of the Model 300 WSL.

A second test was performed using a standard activated charcoal filter. The air flow was measured to be approximately 187.5 cfm. The experiment was set up using the same parameters as outlined above. All samples, including a blank, were sent to an AIHA-certified laboratory for analysis using HPLC and UV detection (NIOSH Method 2016).

When the results of the first test were reviewed, it was discovered that the inlet concentration of formaldehyde during the first test was an order of magnitude less than was found during the second test. The cause of this difference is unknown, but the difference in airflow is believed to have been a factor. Because of the low concentration found during the first test, it was decided to run the test again using an air flow of 187.5 cfm. Accordingly, on April 27, 2011, this third test was performed using an activated charcoal filter that was specially treated for the removal of aldehydes. This time, 2.5 ml of formalin was evaporated into the air stream over a 20-minute period. Air samples were collected at the inlet and outlet of the machine using a sampling rate of about 150cc/min. In addition, an area sample was taken of the room air, placed approximately six feet away from the operation.

Results: The sampling results are presented in the following tables:

**Table I**

**Analysis of Samples collected for Formaldehyde while using the Model 300 WSL  
April 8, 2011; (Using Specially treated Charcoal)**

<b>Sample #</b>	<b>Sample Description</b>	<b>Avg. Flow Rate, cc/min</b>	<b>Sampling time, min.</b>	<b>Sample vol., liters.</b>	<b>Formaldehyde found, mg (ppm)</b>	<b>% Removal</b>
RFA-001	<i>Winged Sentry™</i> Inlet	105.03	30	3.151	0.0025 (0.65)	-----
RFA-002	<i>Winged Sentry™</i> Outlet	104.96	30	3.149	<0.0002 (<0.06)	>96.0

**Table 2**

**Analysis of Samples collected for formaldehyde while using the Model 300 WSL  
April 8, 2011; (Using Standard Activated Charcoal)**

<b>Sample #</b>	<b>Sample Description</b>	<b>Avg. Flow Rate, cc/min</b>	<b>Sampling time, min.</b>	<b>Sample vol., liters.</b>	<b>Formaldehyde found, mg (ppm)</b>	<b>% Removal</b>
RFA-003	<i>Winged Sentry™</i> Inlet	108.98	30	3.151	0.022 (5.5)	-----
RFA-004	<i>Winged Sentry™</i> Outlet	109.32	30	3.149	0.0010 (0.65)	95.45
RFA-005	Blank	----	----	----	<0.0002	-----

**Table 3**

**Analysis of Samples collected for Formaldehyde while using the Model 300 WSL**

**April 27, 2011; (Using Specially treated Charcoal)**

<b>Sample #</b>	<b>Sample Description</b>	<b>Avg. Flow Rate, cc/min</b>	<b>Sampling time, min.</b>	<b>Sample vol., liters.</b>	<b>Formaldehyde found, mg (ppm)</b>	<b>% Removal</b>
RFA-001	<i>Winged Sentry™</i> Inlet	147.04	20	2.941	0.015 (4.2)	-----
RFA-002	<i>Winged Sentry™</i> Outlet	159.26	20	3.185	<0.0002 (<0.06)	>99.33
RFA-003	Area Room Air	154.40	20	3.088	<0.0002 (<0.06)	-----
RFA-004	Blank	-----	-----	-----	<0.0002	-----

**Results and Conclusions:**

1. Both the specially treated activated carbon and standard activated carbon filters used in the Sentry Air Systems SS-300-WSL unit did an excellent job of removing formaldehyde from the air in the experiments described herein. The calculated efficiency of removal ranged between 95.45% for the standard activated carbon to >99.33% for the treated activated carbon.
2. The concentration of formaldehyde in the ambient air was maintained below its detection limits. The Model 300-WSL provided excellent protection throughout the experiment(s). Potential exposures, without regard to the use of a respirator, were found to be well below OSHA workplace limits (PEL) for formaldehyde.

## **Results and Conclusions, continued:**

3. It should be noted that there was some formaldehyde gas found in the sampling tube placed at the outlet of the unit when using standard activated carbon. However, in both experiments where the specially treated activated carbon was used, no formaldehyde was detected in the sampling tubes at the outlet. Therefore, Sentry Air Systems recommends the use of the specially treated activated carbon in the Model 300-WSL and in other Sentry Air System extractors in laboratories and other workplaces for the control of formaldehyde gas.

If you have any questions or comments regarding this report, please contact me at 713-983-7910 or by e-mail at BobCIH@aol.com.

Respectfully submitted,

Robert F. Adams, Industrial Hygienist

RF Adams & Associates

**NOTE: An MSDS (Material Safety Data Sheet) for formalin as used in this project can be found at: <http://avogadro.chem.iastate.edu/MSDS/formaldehyde.htm>**